hydrosense

Hydrosense Ultra Legionella Swab Kit

Instruction Leaflet

SKU 100144

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
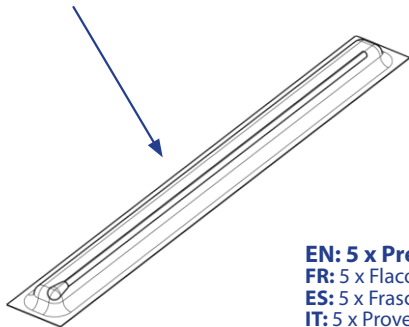
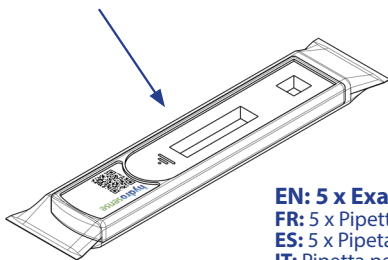
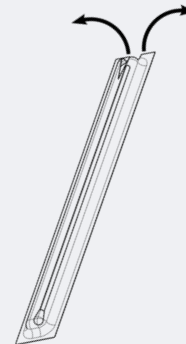
2

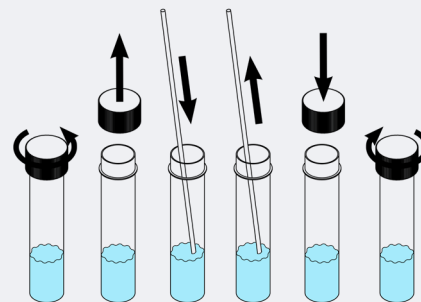EN: If testing a dry surface, pre-moisten the swab by dipping it into the recovery buffer vial.

FR: Si vous testez une surface sèche, pré-humidifiez l'écouvillon en le plongeant dans le flacon de solution de réactif.
ES: Si va a probar una superficie seca, humedezca previamente el hisopo sumergiéndolo en la disolución del frasco.
IT: Se il test viene eseguito su una superficie asciutta, inumidire innanzitutto il tampone immergendolo nella provetta contenente la soluzione tampone.
DE: Wenn Sie eine trockene Oberfläche testen, befeuchten Sie den Tupfer durch Eintauchen in die Wiederherstellungspufferflasche.

3

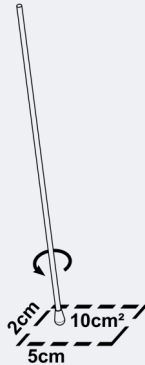
EN: Swab a minimum area of 10cm². Rotate the swab while swabbing the area to achieve more coverage.

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
4

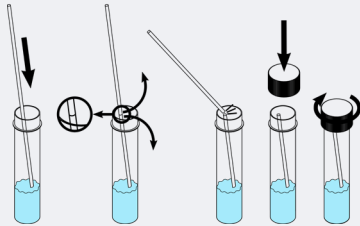EN: Insert swab into the recovery buffer vial. Snap off the swab handle carefully and replace lid.

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5

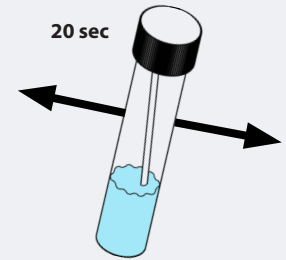
EN: Shake the recovery buffer vial for a minimum of 20 seconds.

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6

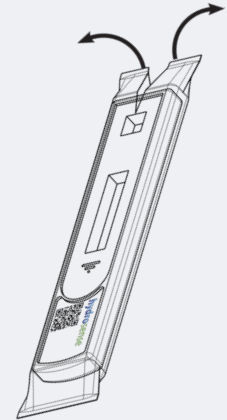
EN: Take the test strip out of its foil packaging. Place the test on a flat surface.

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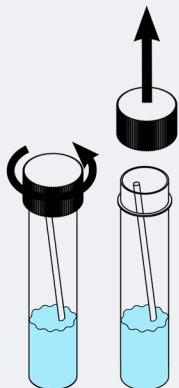
8

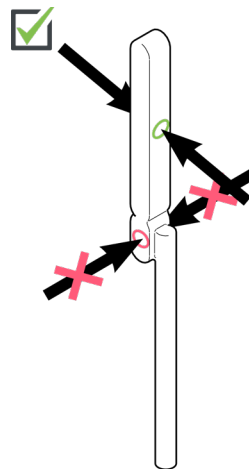
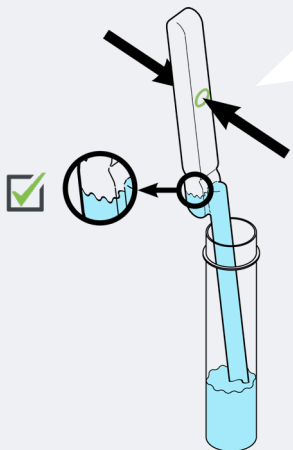
EN: Take exact volume pipette and place tip into recovery buffer solution.

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EN: Press the top bulb ONLY to draw up 0.1 ml of sample into the pipette. Any excess liquid collected in the second (lower) bulb can be ignored.

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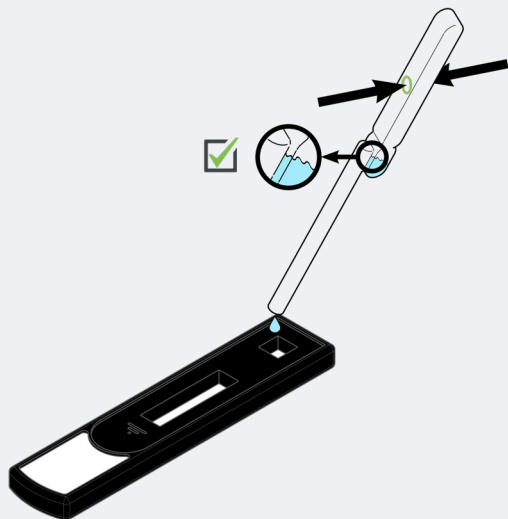
9

EN: Press the top bulb on the pipette to release 0.1ml of sample onto the sample window of the test strip.

Ignore excess in the second (lower) bulb.

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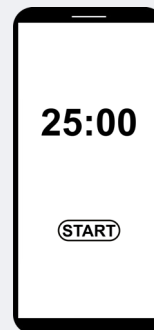
11

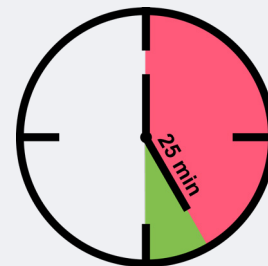
EN: After 25 minutes (and before 30 minutes) read your Hydrosense Legionella test.

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12

EN: The test is composed of two lines; a test line (TL) and a control line (CL). If both lines are present, even if the test line is very faint, then *Legionella pneumophila* serogroup 1 has been detected above the cut-off of the test kit (>200 CFU/ per swabbed area).

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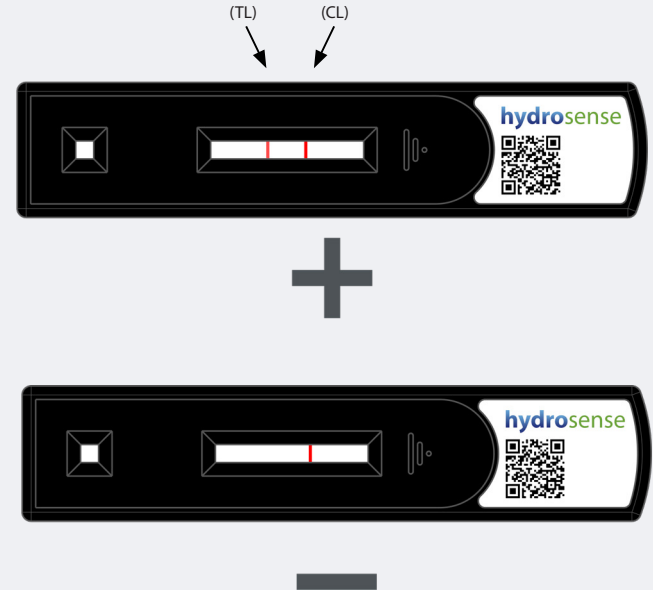
EN: If no test line is present at the test line position but there is a control line, then the test is negative and *Legionella pneumophila* serogroup 1 has not been detected.

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13

EN: Tests without a control line, or that look like any of these tests, are invalid.

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14

EN: Launch the App (Available from Google Play and App Store).

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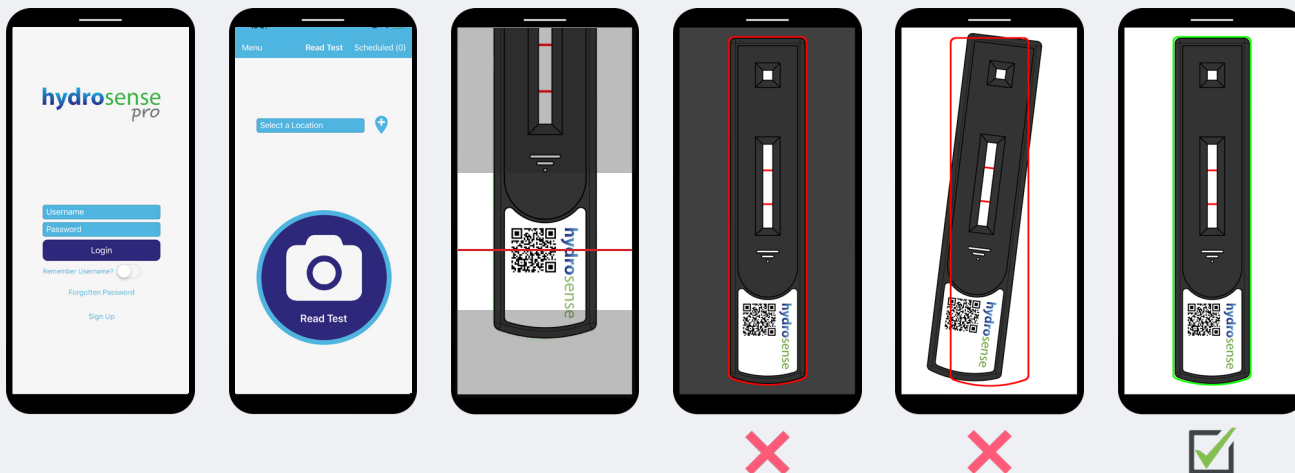
EN: Select a pre-programmed location and click read test. Scan the barcode. The scanner will appear on screen. Read the test by matching up the outline of the test Hydrosense on the screen with the cassette.

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EN: The App will provide you with a positive or negative result and a Hazard index from 0-10. Results can be shared via email and SMS. Results are automatically uploaded to the Hydrosense Pro portal at www.hydrosensepro.com.

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200 KBE *L. pneumophila* sg1 pro Testfläche.

EN: Test Operating Limits:

- Use on samples between 10-45°C (50-113°F).
- Do not use on systems treated with biguanide based or THPS biocides.

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EN: Specificity:

- Non-reactive with wide range of environmental bacteria
- Weak cross reactions with other *L. pneumophila* sg's and *S. aureus* at high concentrations.

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EN: Storage:

- Store at room temperature (18-22°C).
- Do not freeze.
- Do not use after expiry date.
- Do not use if test wrapper perforated.

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IT: Smaltimento

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EN: Disclaimer:

Hydrosense Ltd makes no warranties or representations regarding performance of the products, or that the products are merchantable or fit for a particular purpose. Hydrosense expressly disclaims all other warranties and representations, express or implied, or which arise by operation of law or otherwise.

HYDROSENSE is a trademark of Hydrosense Limited, used worldwide and registered in various territories.

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[©Hydrosense Ltd](http://Hydrosense Ltd)
HYDROSENSE es una marca comercial de Hydrosense Limited, utilizada en todo el mundo y registrada en varios territorios.
[©Hydrosense Ltd](http://www.hydrosense.com)