

Tensormeter RTM1 Product Overview




- ▶ Tensormeter device front panel with 8 signal connectors, reference and trigger connector

- ▶ Simultaneously determine Sheet & Hall resistance at highest precision and extremely low noise
- ▶ Measure irregularly shaped samples without need for lithographic patterning
- ▶ Replace several other devices (Lock-in Amplifier, Source-Measure-Unit, Digital Multimeter, Analog Matrix Switch)
- ▶ Save measuring time, achieve higher throughput




Materials Research and Characterization

- ▶ High precision to study small effects
- ▶ Flexibility for custom measurement sequences
- ▶ Controlled sourcing
- ▶ 2D materials
- ▶ Magnetic materials
- ▶ Transverse resistance

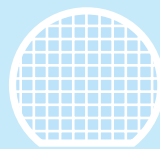

Improved Wafer and Device Testing


- ▶ High stability
- ▶ Faster binning
- ▶ Tighter specs
- ▶ DC and AC
- ▶ Fewer contacts
- ▶ Integrated calibration

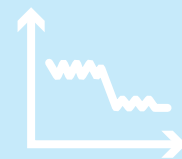

Unique Measurements

- ▶ Irregular sample shapes
- ▶ van-der-Pauw
- ▶ Zero-offset Hall
- ▶ Dummy compensation
- ▶ Device differential
- ▶ Pulse and measure
- ▶ Custom protocols
- ▶ Up to 8 contacts

Interface

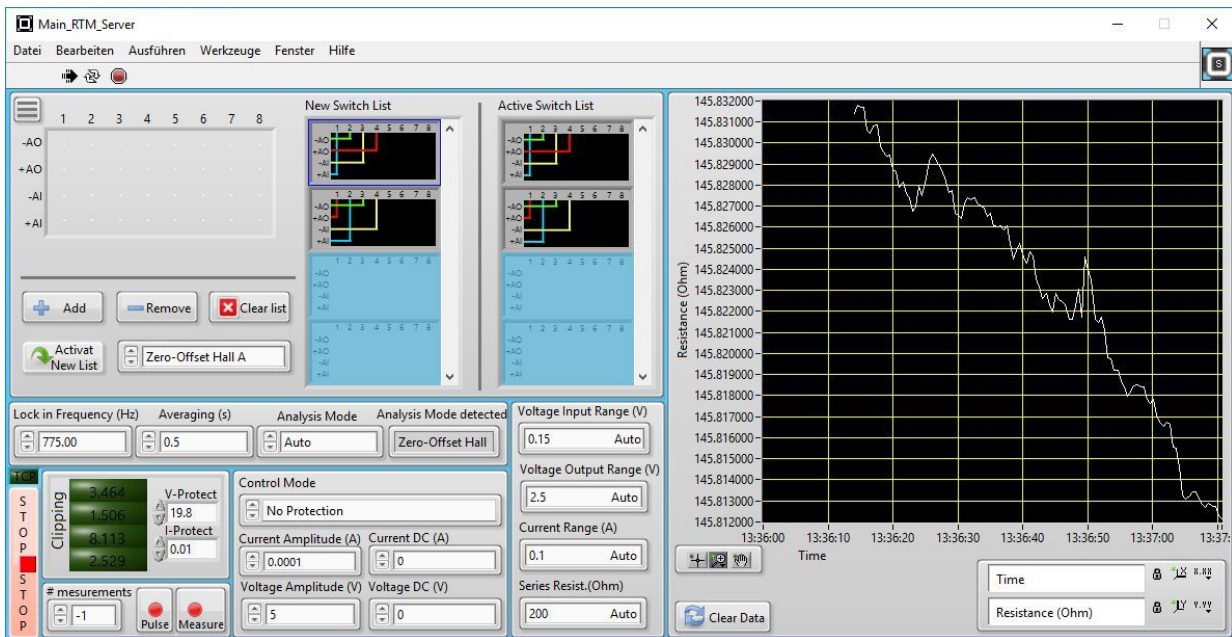

- ▶ Graphical user interface of the server background program, which relays communications between the Tensormeter and the user.

Electrical Specifications

Sensing precision:	<0.1 ppm
Continuous dynamic range:	>8 digits
Symmetrical output:	DC – 20 kHz, ±20 V, ±100 mA
Output noise floor:	< -140 dBFS
Pulse and arbitrary function output with 10 µs resolution	
Input demodulation at multiple frequencies up to 20 kHz	
Differential input noise:	3 nV/√Hz 500 fA /√Hz
Differential input bias current:	1 nA
Optional input transformer for sub-nV/√Hz measurements on low-R DUTs	
Gain change with temperature:	100 ppm/K <1 ppm/K (ratiometric)
DC offset voltage change with temperature:	1 µV/K
Fully controllable integrated 8x4 switching matrix	
Arbitrary function reference input/output:	single-ended ±10 V
Trigger input/output:	single-ended 5V TTL

Software and Communication Protocol Specifications

TCP-based user connection independent of platform and software
Client communication examples for LabView and Python (more on request)
Tensormeter RTM1 connects via USB2.0 to a Windows-PC Software and drivers are provided as Windows Executable Installer
For target OS other than Windows, a small relay computer can be provided
All functions can be controlled from the GUI or via TCP

Hardware, Power and Environmental Specifications


19" rack-mountable device, 3 height units, 25 cm depth
Power demand < 30 W, PSU included, user-specified AC connector
Operation range: 0 – 70 °C, non-condensing humidity
Free convection cooling (can be closed at expense of warmup time)
All front connectors are BNC, 50 Ω type
USB Type B communication connector
Channel and power LED indicators are user-dimmable or can be switched off

Typical Measurement Examples

- ▶ Low noise AC & DC 4-wire measurements in standard geometries (Kelvin, Hall layouts)
- ▶ Presets for van-der-Pauw switched connection 4-wire measurements
- ▶ New Zero-Offset Hall 4-wire preset grants independent longitudinal and transverse resistance
- ▶ Ultra-low noise and high stability Hall measurements outclassing other equipment
- ▶ Sub-ppm relative resistance change investigations
- ▶ Eliminate sample & device drifts with ratiometric resistance measurements
- ▶ High drive harmonic distortion measurements, Pulse & Measure routines, Custom presets



Low Resistive Sensors and Specimen

Differential Input Noise Spectrum of a resistive sensor. Ultra-low wideband & 1/f noise AC measurements allow accurate sensor characterization and operation.

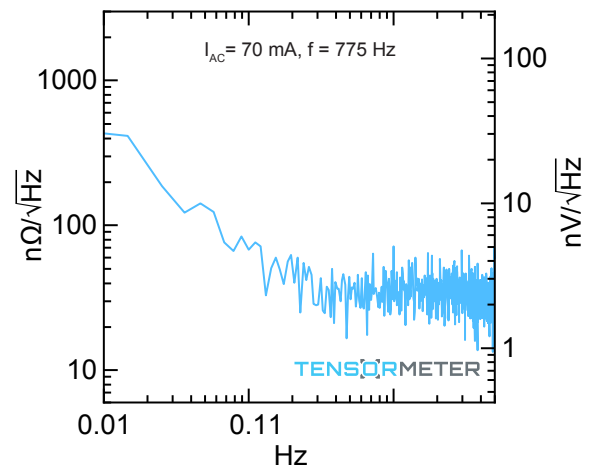


Zero-Offset Hall: Eliminate Drift and Parasitics

Differential Input Noise Spectrum of a Hall measurement on a thin film sample. The Zero-Offset Hall preset of the RTM1 eliminates thermal drift and allows long integration and orders of magnitude improved sensitivity compared to regular 4-wire Hall measurements.

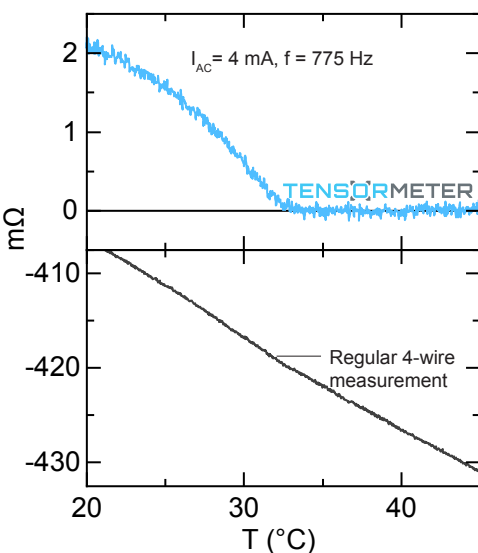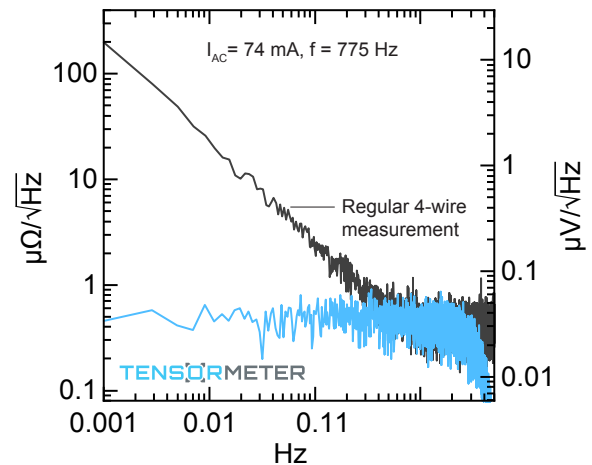

- ◀ Loss of magnetization during warmup of an anti-ferromagnetic sample monitored in Hall Resistance. The Zero-Offset Hall preset of the RTM1 (top) clearly shows the loss of signal. On the contrary, parasitic signal contributions overshadow the useful magnetization signal in a regular 4-wire Hall measurement of the same sample (bottom).

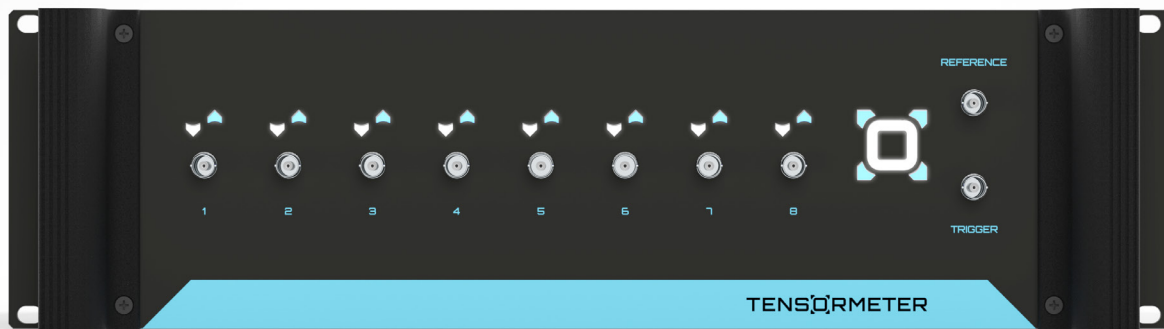


Our Offer



Test

Send us your sample for first test measurements

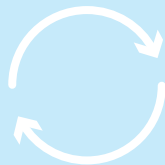

Rent

Try Tensormeter in your lab for some weeks



Buy

Get unlimited Tensormeter performance

Get in Touch

www.tensormeter.de
tensometer@hzdri.de

HZDR Innovation GmbH
Bautzner Landstraße 400
01328 Dresden, Germany

Tensormeter is subject of patents by
Helmholtz-Zentrum Dresden-Rossendorf e.V

HZDR
HELMHOLTZ ZENTRUM
DRESDEN ROSSENDORF

2020 - 03 - 04